- Transition guide/coaxial
- Waveguide-to-coax transition element
- Hohlleiter-Koax-Übergang
- Transizione guida/coassiale
- Transición guía/coaxial

ORITEL TGN 100

FRANÇAIS
ENGLISH
DEUTSCH
ITALIANO
ESPAÑOL

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Meaning of the symbol

CAUTION! Consult the operating manual before using the component.

In this operating manual, the instructions preceded by this symbol may cause personal injury or damage the component and installations if they are not scrupulously abided by or carried out.

You have just bought an **ORITEL TGN 100 waveguide-to-coax transition element** and we thank you for the interest you show for our products.

In order to get the best use out of your component:

- carefully **read** this operating manual,
- **follow** the precautions for use.

PRECAUTIONS FOR USE

- For the waveguide to be terminated correctly, the «N»-type connector of the waveguide-to-coax transition element must always be loaded by its characteristic impedance.
- This component's mass shall be earthed.
- All the instruments and components which are connected to the bench and the AC network shall be earthed.
- No voltage of more than 30 V in relation to the earth shall be present on the bench.

TO ORDER

ORITEL TGN 100 waveguide-to-coax transition element P01.**2753.14**
Delivered with this operating manual

Accessories

ORITEL DEN 100 coaxial detector P01.**2753.15**

CONTENTS

1. PRESENTATION	7
2. DESCRIPTION	7
3. USE	7
3.1 Summary list of the components making up the ORITEL BDH R100 basic educational bench	7
3.2 How to use the EASYFIX quick fastening lug	8
4. CHARACTERISTICS	8
5. MAINTENANCE	9
5.1 Cleaning	9
5.2 Metrological inspection	9
6. GUARANTEE	9
7. ANNEX	22
7.1 General view of the waveguide-to-coax transition element	22
7.2 EASYFIX quick fastening lug	22

1. PRESENTATION

The **ORITEL TGN 100 waveguide-to-coax transition element** is a microwave component specially designed for the ORITEL BDH R100 educational bench, manufactured using a rectangular waveguide in accordance with the R100/WR90 standard (covering the frequency range between 8.5 and 9.6 GHz). This element enables a transmission line in a waveguide to be connected to a female N-type coax and vice versa. By connecting an ORITEL DEN 100 coaxial detector to its output, you have an assembly making it possible to convert the modulated microwave signal to weak or zero frequency signals.

2. DESCRIPTION

See § 7.1

- ① Female «N» connector.
- ② Knurled tightening knob for the EASYFIX quick fastening lug.

3. USE

Generally speaking, the components below are a part of the standard supplies delivered with the ORITEL BDH R100 basic educational bench.
To use these components, refer to their respective operating manuals.

3.1 Summary list of the components making up the ORITEL BDH R100 basic educational bench

ORITEL OSG 100 Gunn diode oscillator	P01.2753.07
ORITEL ISO 100 ferrite isolator	P01.2753.08
ORITEL MOD 100 pin diode modulator	P01.2753.09
ORITEL ATM 100 micrometer-adjustable variable attenuator	P01.2753.10
ORITEL OND 100 curve wavemeter	P01.2753.11
ORITEL LAF 100 measuring line	P01.2753.12
ORITEL ADZ 100/3 three-screw impedance adaptor	P01.2753.13
ORITEL TGN 100 waveguide-to-coax transition element	P01.2753.14
ORITEL DEN 100 coaxial detector	P01.2753.15
ORITEL CHG 100 adapted load	P01.2753.16
ORITEL CC 100 short circuit plate	P01.2753.17
ORITEL SUP 100 waveguide support	P01.2753.18

■ Other components (R100/WR90 waveguide)

ORITEL AFR 100 EASYFIX quick fastening lug adaptor	P01.2753.01
ORITEL RS 100 bench support rail	P01.2753.03
ORITEL MH 600 microwave milliwattmeter	P01.2501.01
ORITEL ST 613 thermocouple probe (for ORITEL MH 600)	P01.2851.01
ORITEL CF 204A power supply unit for Gunn diode oscillator	Please consult us
ORITEL IR 205 SWR indicator	P01.2705.01
ORITEL RD 100 displacement sensor for measuring line	P01.2753.02
ORITEL ANC 100/15 (15 dB) horn antenna	P01.2753.04
ORITEL CGX 100/20 waveguide cross coupler	P01.2753.05
ORITEL IRIS 20-30 (20 and 30 dB) coupling irises	P01.2753.06

To use these components, refer to the corresponding operating manual.

3.2 How to use the EASYFIX quick fastening lug

See diagram § 7.2

- Arrange the two elements to be assembled, taking care to correctly place the centring studs ① opposite the holes ②
- Simultaneously tighten the two knurled knobs ③

4. CHARACTERISTICS

Frequency: 8.5 to 9.6 GHz

SWR: < 1.25 of 8.5 to 9.6 GHz

Waveguide: R100

Flange: UBR 100

Connector: female «N»

Weight: 180 g

Dimensions: 56.5 x 53 x 41.5 mm

Electromagnetic compatibility:

Compliance with the NF EN 61326-1 standard, 1998 edition

5. MAINTENANCE

⚠ **For maintenance, use only specified spare parts. The manufacturer will not be held responsible for any accident occurring following a repair done other than by its After Sales Service or approved repairers.**

5.1 Cleaning

Before assembling the educational bench's microwave components, make sure that no dust is present inside the waveguide.

⚠ **CAUTION!**
Do not use a pressurised air jet for fear of damaging certain components.
Do not use solvent, and take precautions when dusting.

5.2 Metrological inspection

⚠ **It is essential that all measuring instruments are regularly calibrated.**
For checking and calibration of your instrument, please contact our accredited laboratories (list on request) or the Chauvin Arnoux subsidiary or Agent in your country.

■ **Repairs under or out of guarantee**
Please return the product to your distributor.

6. GUARANTEE

Our guarantee is applicable for **twelve months** after the date on which the equipment is made available (extract from our General Conditions of Sale, available on request).

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